

Stepper StepMate

Features:

- ◆ Light and ergonomic solution provides easy usability
- ◆ Single-handed operation
- ◆ Volume range from 10µL to 5mL
- ◆ Up to 48 dispensing steps
- ◆ Maintenance free
- ◆ Equipped with durable tip

Optimized for smooth dispensing of even the smallest volumes

Five easy-to-select settings control dispensing volume

How to select dispensing volume

Find the required dispensing volume in the table
Use the adjusting wheel to set the required volume
Select and insert a suitable tip
Follow the maximum dispensing steps

Tip volume		Dispensing Volume				
Tips	Sel.	1	2	3	4	5
0.50ml		10µl	20µl	30µl	40µl	50µl
1.25ml		25µl	50µl	75µl	100µl	125µl
2.50ml		50µl	100µl	150µl	200µl	250µl
5.00ml		100µl	200µl	300µl	400µl	500µl
12.50ml		250µl	500µl	750µl	1000µl	1250µl
25.00ml		500µl	1000µl	1500µl	2000µl	2500µl
50.00ml		1000µl	2000µl	3000µl	4000µl	5000µl
Steps		48	23	15	11	8



By pushing the slide upwards, the syringe is filled up

The stop lever secures the syringe

Dial setting		1	2	3	4	5			
Number of steps		48	23	15	11	8			
Cat. No.	Syringe volume (mL)	Sample volume (µL)					A %	P %	Description
17900032	0.50	10	20	30	40	50	± 0.8	≤ 0.7	0.5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900033	1.25	25	50	75	100	125	± 0.8	≤ 0.5	1.25mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900034	2.50	50	100	150	200	250	± 0.8	≤ 0.4	2.5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900035	5.00	100	200	300	400	500	± 0.4	≤ 0.3	5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900036	12.50	250	500	750	1000	1250	± 0.3	≤ 0.3	12.5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing



Dial setting		1	2	3	4	5			
17900037	25.00	500	1000	1500	2000	2500	± 0.2	≤ 0.3	25mL, piston PE-HD, cylinder PP, 1 packing, 2pcs/packing with 1 pc adapter
17900038	50.00	1000	2000	3000	4000	5000	± 0.2	≤ 0.2	50mL, piston PE-HD, cylinder PP, 1 packing, 2pcs/packing with 1 pc adapter